



Syrris Ltd.
27 Jarman Way, Royston, Hertfordshire
SG8 5HW, United Kingdom
T: +44 (0)1763 242555
F: +44 (0)1763 242992
E: info@syrris.com
W: www.syrris.com

Syrris Limited
Registered office as address
Company No. 4257809

| | |
|------------------------|---|
| System/Product | Atlas XL Pump Valve Syringe Compatibility |
| Model/Module | Atlas |
| Subject Type | Sales training/update to instruction manual |
| Subject Version | |
| Author | RJR |
| Date | 14 th December 2009 |

The following table highlights the compatibility between syringes and valves for the Atlas XL pump.

Key

- ✓ No Conflict
- X Port positions blocked (no suitable fittings)
- X¹ Port positions partially blocked (cannot use large head gripper, can use compact gripper and ferrule type)
- X² Port positions partially blocked (cannot use large head gripper or ferrule type, can use compact gripper)

| | 3-Way | 4-Way | 6-Way | 8-Way |
|--------------|--------------|--------------|----------------|----------------|
| 500ul | ✓ | ✓ | D ¹ | E ¹ |
| 1ml | ✓ | ✓ | D ¹ | E ¹ |
| 2.5ml | ✓ | ✓ | D ² | E ² |
| 5ml | ✓ | ✓ | D | E |
| 10ml | ✓ | ✓ | D | E |
| 25ml | ✓ | ✓ | D | E |
| 50ml | ✓ | C D | D | D E F |